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• **Sample preparation:**

**High quality DNA is essential for high quality results** (For template re-purification an extra cost of 5.00 Euro (SADN 0004) will be charge). We recommend to use mini or midi prep with ion exchange resin purification (Quiagen, Nucléospin, GibcoBRL Concert, Sigma,...). Dilute DNA **in water** (avoid TE buffer) and check the concentration with a spectrophotometer and DNA integrity by electrophoresis of 3µl on agarose gel and **please send the picture**. Send 2-3µg of plasmid (**0.2 µg/µl**) in 15 µl of **pure water (avoid TE)**. If your DNA contains a special specificities (high GC content, repeats, homopolymic stretches,...) please order SADN 0007.

• **PCR products:**

Send the purified (20ng/100 bases and **resuspend in pure water, avoid TE or any buffer**) or non-purified PCR product (order SPCR 0002 or SPCR0005 if several products) and your primer(s) (10 µM) or the primer sequence(s) (order SSOL 0003).

• **Confidentiality:**

We guarantee to maintain the confidentiality of our data at any time. 3 months after delivery your complete data will be deleted and residual DNA will be dismissed.

• **Order:**

Please use the DNA sequencing order form that you can download from our web site (<http://www.millegen.com>) and send it with the purchase order of your institute (if you cannot send them simultaneously, please **indicate the purchase order number (PO #)** on the DNA sequencing order form) and the sample to our address:

**MilleGen**, Prologue Biotech, rue Pierre et Marie Curie, BP 700, 31319 Labège (France) Tel:+ 33 (0)5 61 28 70 16 ; Fax: 33 (0) 5 61 28 70 11; Email: [sequence@millegen.com](mailto:sequence@millegen.com)